




Hi-Temp Oven

Model: DR210C

- First edition -

- Thank you very much for purchasing this Yamato DR210C Hi-Temp Oven.
- Please read the “Operating Instructions” and “Warranty” before operating this unit to assure proper operation. After reading these documents, be sure to store them securely together with the “Warranty” at a handy place for future reference.
-  **WARNING!:**
Carefully read and thoroughly understand the important warning items described in this manual before using this unit.

Yamato Scientific Co.,Ltd.

Table of contents

1. Safety precautions	1
Explanation of pictograms	1
List of symbols.....	2
Warning • Cautions	3
2. Before operating the unit	4
Precautions when installing the unit	4
3. Names and functions of parts	8
Main body	8
Operation panel.....	9
Explanation of characters	10
4. Operating procedures	12
List of operation modes and functions.....	12
Operation mode • function setting keys and characters	14
Settings for overheat prevention device	15
Operating procedures (fixed temperature operation).....	16
Operating procedures (quick auto stop operation)	17
Operating procedures (auto stop operation).....	19
Operating procedures (auto start operation).....	21
Preparing a program.....	23
Program repeat operation.....	28
Programming sheet.....	29
Useful functions (calibration offset function)	31
Useful function (lock function)	32
5. Cautions on handling	33
6. Maintenance procedures	36
Daily inspection/maintenance	36
7. When the unit is not to be used for a long time or when disposing	37
When the unit is not to be used for a long time or when disposing.....	37
Notes about disposition	37
8. Troubleshooting	38
Safety device and error codes.....	38
When a malfunction is suspected	39
9. After sales service and warranty	40
When requesting a repair	40
10. Specifications	41
11. Wiring diagram.....	42
12. List of replacement parts	43
13. List of dangerous materials.....	44
14. Standard installation manual	45

1. Safety precautions

Explanation of pictograms

About pictograms

A variety of pictograms are indicated in this operating instruction and on products to assure safe operation. Possible results from improper operation ignoring them are classified as follows.

Be sure to fully understand the descriptions below before proceeding to the text.



Warning

Indicates a situation which may result in death or serious injury (Note 1)



Caution

Indicates a situation which may result in minor injury (Note 2) and property damage (Note 3).

(Note 1) Serious injury means a wound, an electrical shock, a bone fracture or intoxication that may leave after effects or require hospitalization or outpatient visits for a long time.

(Note 2) Minor injury means a wound or an electrical shock that does not require hospitalization or outpatient visits for a long time.

(Note 3) Property damage means damage to facilities, devices and buildings or other properties.

Meanings of pictograms



This pictogram indicates a matter that encourages the user to adhere to warning ("caution" included).
Specific description of warning is indicated near this pictogram.



This pictogram indicates prohibitions
Specific prohibition is indicated near this pictogram.



This pictogram indicates matters that the user must perform.
Specific instruction is indicated near this pictogram.


1. Safety precautions

List of symbols

Warning



General warnings

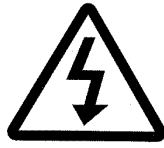

Danger!:
High voltage



Danger!:
High temperature



Danger!:
Moving part



Danger! Hazard of
explosion

Caution



General cautions



Electrical shock!



Burning!



Caution for no
liquid heating!



Caution for water
leak!



For water only



Poisonous
material

Prohibitions



General bans



Fire ban



Do not
disassemble



Do not touch

Compulsions



General
compulsions



Connect ground
wire



Install levelly



Pull out the power
plug



Regular
inspection

1. Safety precautions

Warning • Cautions

Warning



Never operate the unit in an atmosphere containing flammable or explosive gas

Never operate the unit in an atmosphere containing flammable or explosive gas. Otherwise, an explosion or a fire may result since the unit is not explosion-proof. See section “13. List of dangerous materials” on page 44.



Be sure to connect the ground wire.

Be sure to connect the ground wire correctly. Otherwise, electrical leak may result and cause an electrical shock or a fire.



Ban on operation when an abnormality occurs

When a smoke or a unusual odor is seen or sensed, immediately turn the power switch on the main unit off and pull out the power cord (plug) from the power supply. A fire or an electrical shock may result.



Never use electrical power cords bundled.

When these are used bundled, they might overheat causing a fire.



Take care not to damage electrical power cords.

Avoid tightly bend, pull with a strong force or twist to prevent electrical power cords from damaging. A fire or an electrical shock may result.



Never use an explosive or a flammable material with this unit.

Never use an explosive material, a flammable material or a material containing them. An explosion or an electrical shock may result. See section “13. List of dangerous materials” on page 44.



Never try to touch a hot part.

Some parts of the unit are hot during and immediately after operation. Take special care for possible burning.



Never try to disassemble or alter the unit.

Never try to disassemble or alter the unit. A malfunction, a fire or an electrical shock may result.



Caution



When a thunder is heard.

When a thunder is heard, turn the main power off immediately. A malfunction, fire or an electrical shock may result.

2. Before operating the unit

Precautions when installing the unit



Warning

1. Be sure to connect the ground wire.



- DR210C types use a single-phase 220V power source. Please consult your local electrical contractor for power connecting work.
- Be sure to connect the ground wire (green core wire of the power cord) to the ground wire or the ground terminal in order to prevent an electrical shock from electric leakage.
- Never connect the ground wire to gas or water pipes. Otherwise, a fire may result.
- Never connect the ground wire to the ground for telephone wire or lightning conductor. Otherwise, a fire or an electrical shock may result.
- Do not use a branching receptacle, which may generate heat and possible danger.

2. Carefully select an installation site.




Take special care not to install the unit at a place described below:

- On uneven or dirty floor
- Where combustible gas or corrosive gas exists
- Where the ambient temperature is 35°C or more
- Where temperature widely fluctuates
- Where there is excessive dust or humidity
- Where subject to direct sunlight
- Where vibration is severe



Install this unit at a place with spaces shown below.
Do not cover the exhaust ports on the top of the unit.

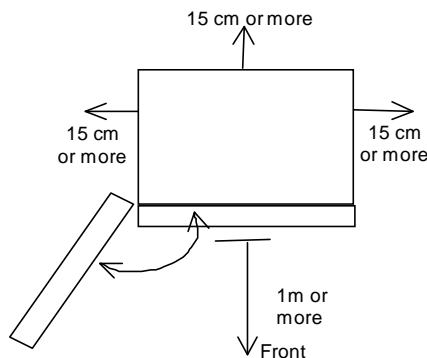




2. Before operating the unit


Precautions when installing the unit


Warning

3. Never operate the unit in an atmosphere containing explosive or flammable gas

-  Never operate the unit in an atmosphere containing flammable or explosive gas. Since the unit is not explosion-proof, an arc is discharged when turning a switch “ON” and “OFF” and during operation and a fire or an explosion may result.
-  See the section “13. List of dangerous materials” on page 44 for flammable and explosive gases.

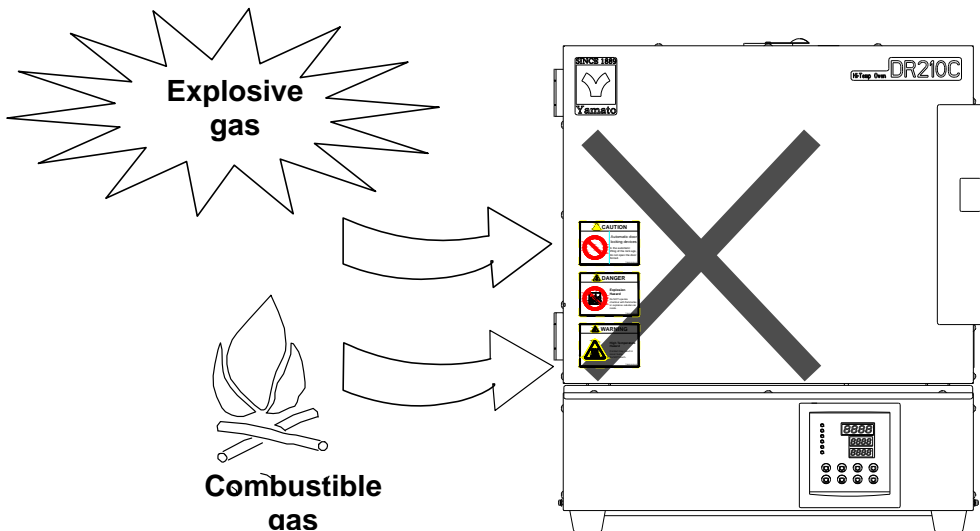

-  Never use the unit with an explosive material, a flammable material or a material containing them. An explosion or a fire may result.
- See the section “13. List of dangerous materials” on page 44 for flammable and explosive substances.

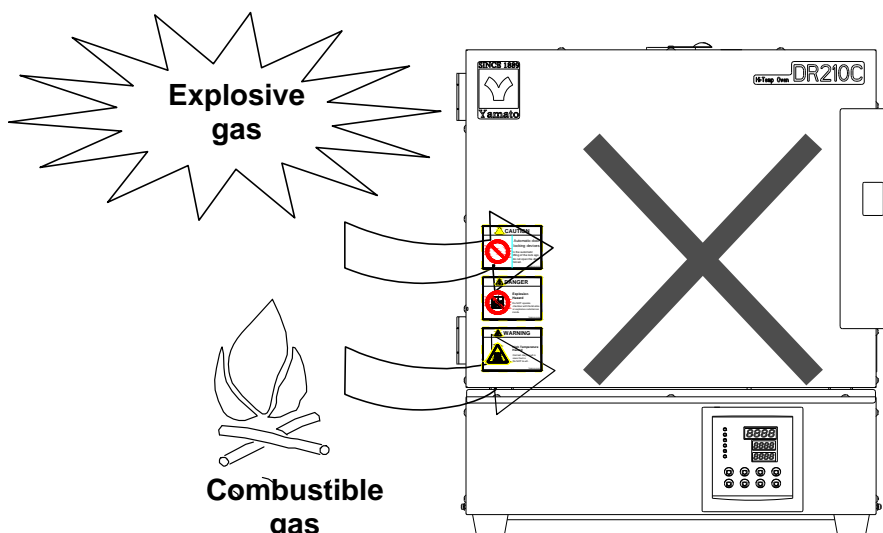


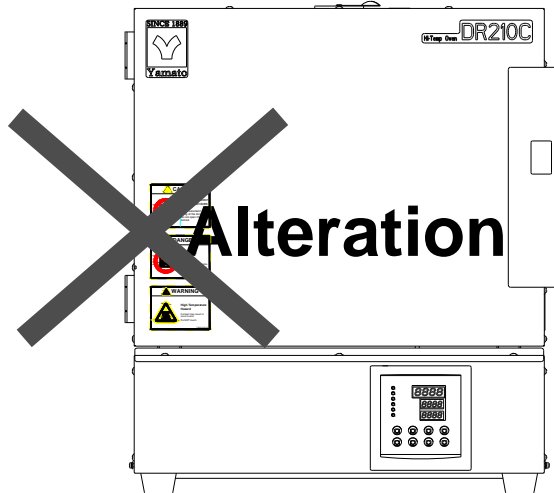
2. Before operating the unit

Precautions when installing the unit


Warning


4. Do not alter the product.

-  The user shall never attempt to alter the unit since it may cause a malfunction.




5. Install the unit on a level surface

-  Install the unit on a level surface. Installing this unit on a slope might cause unexpected troubles or malfunctions.

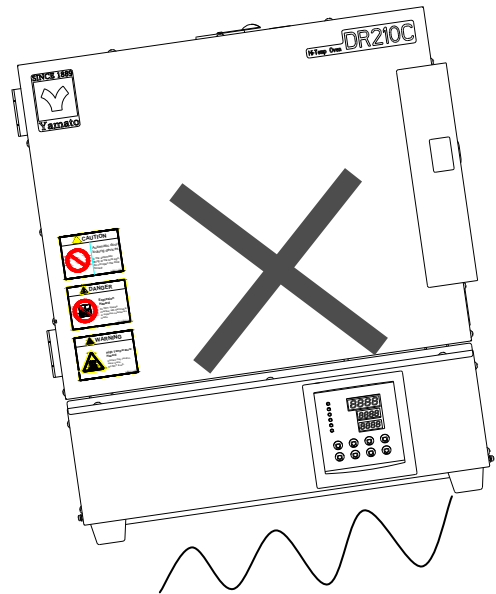


6. Be sure to connect the power plug to the dedicated wall outlet.

-  Use a power distribution panel or a wall outlet that meets the electrical capacity of the unit.

Electric capacity: DR210C AC220V 6A

When the unit will not start even when you turn the power switch to "ON", check for low main voltage or if the unit is connected to the same power supply line as other devices and connect it to another line if necessary.

7. Installation

-  The unit might fall down or move by an earthquake or an impact resulting a personal injury. We recommend making safety measures such as to avoid installing the unit at a place other than busy places.
Also, take appropriate safety measures to prevent the unit from tripping over.

2. Before operating the unit

Precautions when installing the unit



Caution

8. About handling of the power cord



Never use electrical power cords bundled. When these are used bundled, they might overheat causing a fire.

Do not convert, forcibly bend, twist or pull the power cord. Otherwise, a fire or an electrical shock may result.

Do not place the power cord under a desk or a chair, or sand between objects to avoid it from being damaged.



Do not place the power cord close to a stove or other heat generating device. Sheath of the cord may burn and result in a fire or an electrical shock.

If the power cord should be damaged (exposure of core wire or disconnection), immediately turn the main unit off, pull out the power cord (plug) out of the power supply and ask your dealer to replace the cord. Otherwise, a fire or an electrical shock may result.

Connect the power cord to an appropriate wall outlet.


9. When you operate the unit for the first time



When you operate the unit for the first time at a higher temperature, the unit may generate an odor. This is due to decomposed bonding material contained in heat-insulation material and is not a malfunction of the unit. We recommend operating the unit at the highest temperature once before starting its regular operation.


3. Names and functions of parts

Main body

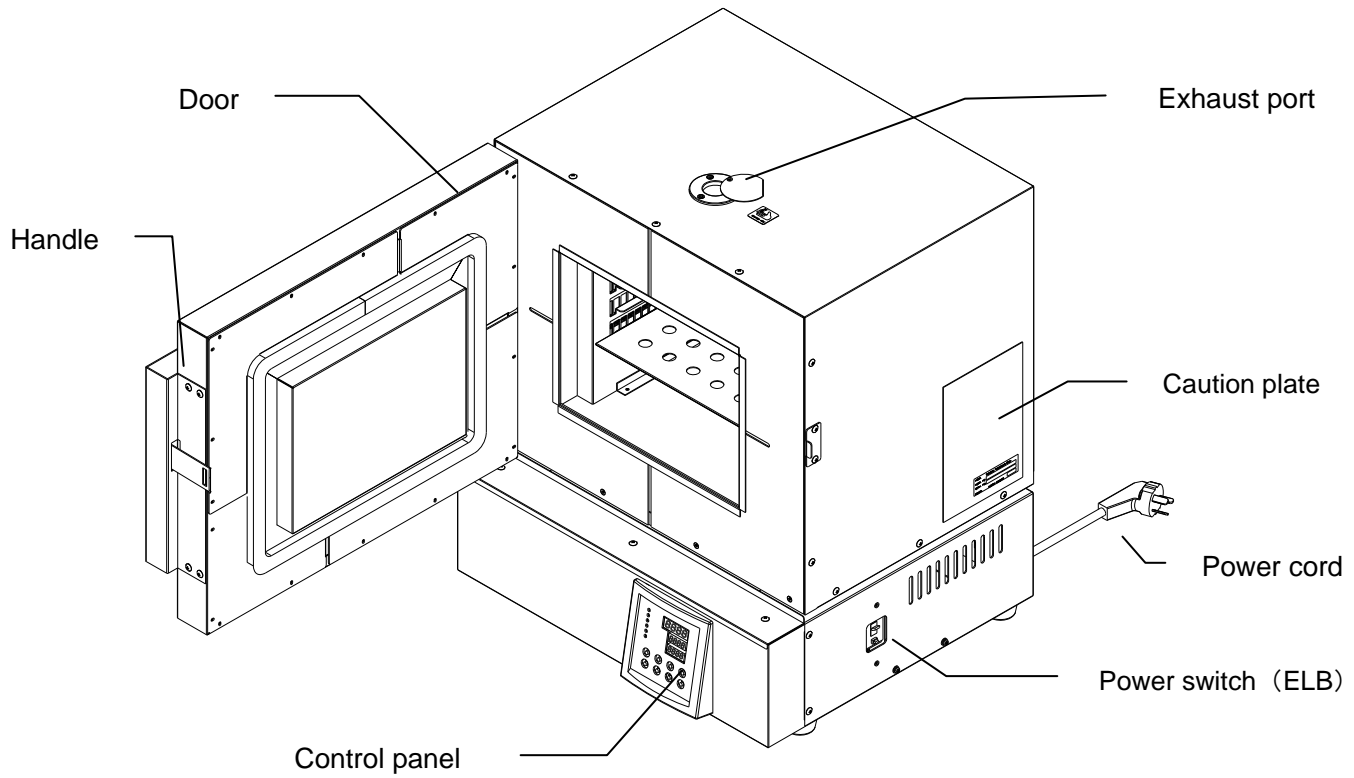

3. Names and functions of parts

Operation panel

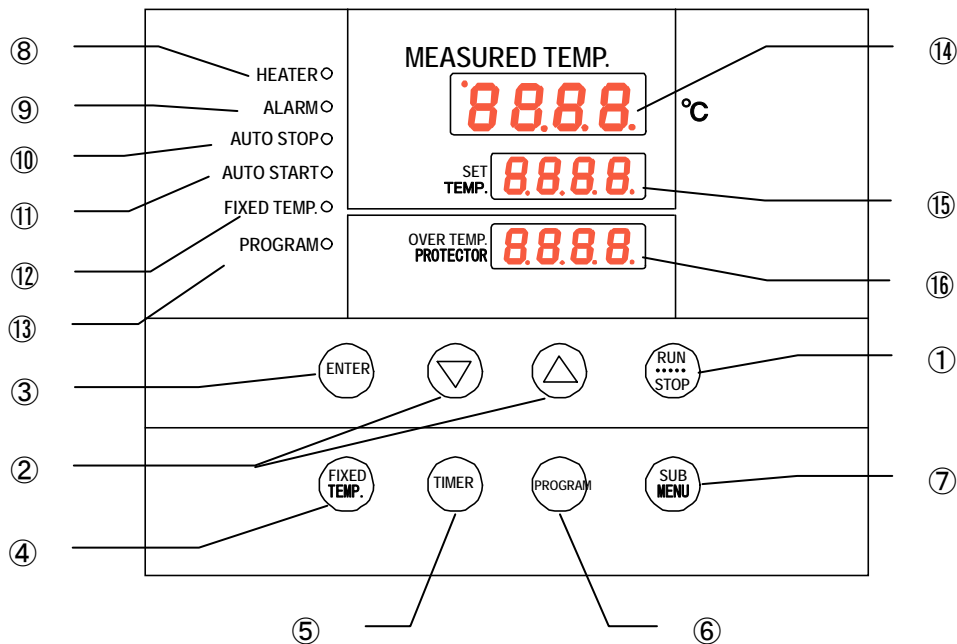






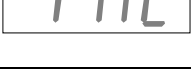



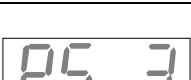


No.	Name	Operation/action
①	RUN/STOP key	Used for starting/stopping operation
②	▲▼ keys	Used for selecting settings.
③	ENTER key	Used for determining a selected setting.
④	FIXED TEMP. key	Key for selecting fixed temperature operation.
⑤	TIMER key	Key for selecting timer operation. Quick auto stop operation, auto stop operation or auto start operation can be selected.
⑥	PROGRAM key	Used for selecting programming and starting the program operation. Programs of six patterns in three types can be set.
⑦	SUB MENU key	Key for setting temperature of the overheat prevention device, calibration offset temperature, the key lock function or the program repeat function.
⑧	HEATER lamp	Illuminates while heater power is on.
⑨	ALARM lamp	When an error occurs, the lamp illuminates with an audible buzzer.
⑩	AUTO STOP lamp	Flashes while the quick auto stop timer is being set and illuminates while the unit is in operation. Flashes while the auto stop timer is being set and illuminates while the unit is in operation.
⑪	AUTO START lamp	Flashes while the auto start timer is being set and illuminates while the unit is in operation.
⑫	FIXED TEMP. lamp	Flashes while fixed temperature operation is being set and illuminates while the unit is in operation.
⑬	PROGRAM lamp	Flashes while the program operation is being set and illuminates while the unit is in operation.
⑭	MEASURED TEMP. display	Displays measured temperature in the bath • set characters • alarm information.
⑮	SET TEMP. display	Displays a set temperature, timer settings and timer remaining time.
⑯	OVER TEMP. PROTECTOR display	Displays the set temperature for the overheat prevention device.

3. Names and functions of parts




Explanation of characters

Characters on the VS6 type controller are explained in this section.

Characters	Identifier	Name	Application
	Fix	Fixed temperature operation setting	Means settings for fixed temperature operation.
	Sv	Temperature setting	Used for setting a temperature.
	AStP	Auto stop setting	Used for setting auto stop operation.
	AStr	Auto start setting	Used for setting auto start operation.
	tim	Time setting	Used for setting a time.
	PrG3	Program type select	Used for selecting a program type 1, 2 or 3 to be used. See section "Preparing a program on page 23."
	PA1	Program pattern select	Used for selecting a program pattern to be used. See section "Preparing a program on page 23."
	End	Time up	Displayed when timer operation has ended. See page 17, 19, and 21.
	Sv-1	Program temperature setting	Used for setting temperatures for each step of a program. (Sv-1 to Sv-30 are displayed.)
	t-1	Program time setting	Used for setting time for each step of a program. (t-1 to t-30 are displayed.)
	PS-3	Program repeat Selecting a return destination	Used for selecting a step for returning during program repeat operation. See section "Program repeat operation" on page 28.
	Pc-2	Setting number of program repetitions	Used for setting number of program repeat operations. See section "Program repeat operation" on page 28.

3. Names and functions of parts

Explanation of characters

Characters	Identifier	Name	Application
	cAL	Calibration offset setting	Used for inputting a calibration offset temperature See section "Using the calibration offset function" on page 31.
	oH	Setting overheat protection device temperature	Used for setting an overheat protection device temperature. See section "Settings for overheat prevention device" on page 15.
	Lock	Key lock of settings	Key locks settings to prevent their alteration See section "Using the lock function" on page 32.

* See the section "Operation mode • function setting keys and characters" on page 14 for characters of operation modes and functions

4. Operating procedures

List of operation modes and functions

Operation modes of the unit are as shown below:

No.	Name	Description	Page
1	Fixed temperature operation	Pressing the FIXED TEMP. key to enter the fixed temperature operation setting mode. Pressing the FIXED TEMP. key again to enter the temperature setting mode. Set a temperature with the ▼▲ keys. Pressing the RUN/STOP key to start operation, and pressing the RUN/STOP key again to stop operation.	P.16
2	Quick auto stop operation	Used when you want to, for example, “stop operation being performed automatically in several hours.” Pressing the TIMER key during the fixed temperature operation to enable setting a time before operation stops. Set a duration before stop with the ▼▲ keys. Pressing the RUN/STOP key starts quick auto stop operation and activates the timer in the middle of it to automatically stop it after the set period of time.	P.17
3	Auto stop operation	Used when you want to “set automatic stop for fixed temperature operation when making settings for it.” Press the TIMER key to display “AStP.” Pressing the ENTER key to set the temperature setting “Sv.” Pressing the ENTER key again to enable setting of the operation time “tim.” Pressing the RUN/STOP key starts auto stop operation.	P.19
4	Auto start operation	Used when you want to “start operation automatically after several hours” after power is turned on. Press the TIMER key to display “AStr.” Pressing the ENTER key to set the temperature setting “Sv.” Pressing the ENTER key again to enable setting of the operation time “tim.” Pressing the RUN/STOP key starts auto start operation.	P.21
5	Program operation	This is used to raise or lower temperature based on the set temperature and time. Press the PROGRAM key to display “PrGn.” (n:1,2,3) Press the PROGRAM key again to select a program mode you want. To select “PrG2” or “PrG3”, press the ENTER key to select a pattern “PAT” you want. Pressing the RUN/STOP key starts program operation.	P.23
* Operation mode cannot be changed while the unit is in operation. First stop operation before changing the mode.			

4. Operating procedures

List of operation modes and functions

Operating functions of the unit are as shown below:

No.	Name	Description	Page
1	Overheat prevention function	<p>Automatic overheat prevention function: This function is linked to the unit set temperature and has been set to so that it is automatically activated (returned automatically) at a temperature 12°C higher than the set temperature in the bath.</p> <p>Overheat prevention device: Although the power supply, the display and the key input assembly are in common with the controller, the device consists of the standalone temperature measurement circuit, the CPU, the sensor, and the output circuit enabling to set to a temperature you want on the operation panel. If the overheat prevention device is triggered, the unit will stop and will not recover until the power switch is turned on again. (Manual recovery)</p>	P.15
2	Calibration offset function	<p>Calibration offset function compensates any differences between the target temperature in the bath and the control temperature of the controller (sensor temperature.) The function can compensate to either plus or minus side for the whole temperature band of the unit.</p>	P.31
3	Overheat prevention temperature compensation function	When temperature compensation in item two is made for the controller, temperature to be input to the overheat prevention device will also be compensated automatically.	—
4	Power outage compensation function	When power outage occurs in the middle of operation, the operation resumes from the status immediately before the power outage.	—
5	Setting lock function	<p>This function locks the set operation status. The lock can be set or released with the SUB MENU key.</p>	P.32

Operation mode · function setting keys and characters

4. Operating procedures

Settings for overheat prevention device

The safety device for preventing overheat has the automatic overheat prevention function for the controller (automatic recovery) and has the power supply, the display and the key input assembly in common with the controller as well as it has an overheat prevention device (manual recovery) as a secondary safety measures that consists of separate temperature measurement circuit, the CPU, the sensor and the output circuit.

Temperature setting range and functions

The device has overheat prevention function in double. One is the function integrated in the controller, which is designed to be activated automatically at a temperature 12°C above the set temperature of the temperature controller (the heater repeats ON/OFF at +12°C) at the time of shipping from the factory.

Another is integrated with the controller and is set with the keys on the controller.


The secondary overheat prevention function is added by this setting.

The temperature range settable for the overheat prevention device integrated with the controller is "0°C to the 50°C + the highest set temperature for the device."

When the temperature in the bath keeps rising beyond the controller set temperature and reaches the set temperature of the overheat prevention device, the circuit trips, Er19 flashes on the controller screen, and the buzzer continues to sounds.

When the overheat prevention device is activated, it will not be released until power is turned on and Er19 will not be released.

How to set temperature

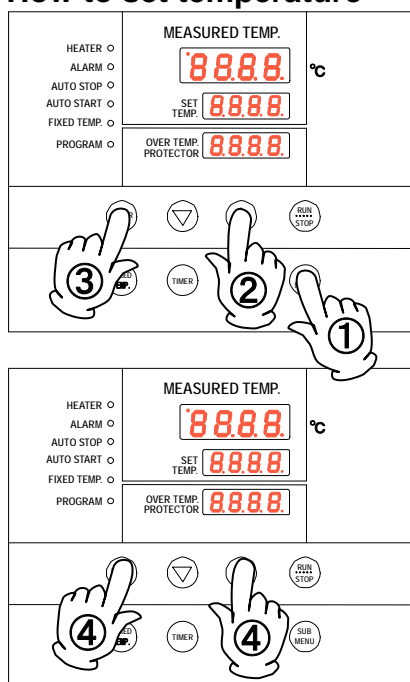

1. Turn power ON. (Turn the ELB to ON.)

When power is turned ON, the initial values will be displayed for about four seconds, then the initial screen will appear and the current bath temperature, operation mode character and the overheat prevention set temperature are displayed on each of the displays.

2. Setting the overheat prevention temperature

- ① Press the **SUB MENU** key.
- ② Press the **▼▲** keys to select the overheat prevention temperature setting character **oH** for the MEASURED TEMP. display.
- ③ Pressing the **ENTER** key displays the current set temperature flashing on the SET TEMP. display.
Caution: In general, set to a temperature at least 10°C higher than the controller set temperature to prevent the device from malfunctioning.
- ④ When the temperature you want is obtained with the **▼▲** keys, press the **ENTER** key to complete setting.


Caution

- ① Set temperature as "highest operation temperature for the unit +30°C" or "set temperature +30°C" as a rough standard and add 5°C to the setting if the device functions improperly.
- ② Be sure to set the overheat prevention activation temperature correctly otherwise the device may not start, the overheat prevention device is activated before temperature in the bath increases completely, or a fire or other unexpected accidents may result.
The temperature is set at 730°C on shipping from the factory.
- ③ The overheat prevention device has been designed to prevent overheating of devices not to protect samples. The device does not prevent accidents caused from use of explosive or flammable materials.

4. Operating procedures

Operating procedures (fixed temperature operation)

How to start fixed temperature operation

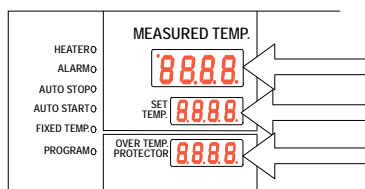

1. Turn power ON. (Turn the ELB to ON.)


When power is turned ON, the initial values will be displayed for about four seconds, then the initial screen will appear and the current bath temperature, operation mode character and the overheat prevention set temperature are displayed on each of the displays.

MEASURED TEMP. display: Indicates the present bath temperature and various characters.

Set temperature display: Indicates the set temperature and various characters.

Overheat prevention set temperature display: Indicates the set temperature for the overheat prevention device.


For operation mode characters, see section page 10.

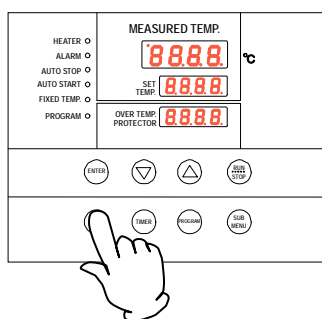

2. Selecting an operation mode

Press the **FIXED TEMP.** key to display the fixed temperature operation mode on the SET TEMP. display.

Display the character Fix **Fi II**.

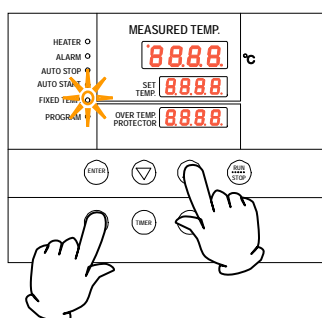


3. Setting the temperature

Press the **FIXED TEMP.** key again.

The character SV **50** that indicate a set temperature is displayed on the MEASURED TEMP. display on which the current set temperature flashes and the FIXED TEMP. lamp flashes.

Set a temperature using the **▼▲** keys.

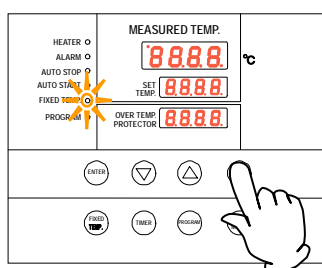

4. Starting operation

Press the **RUN/STOP** key for about one second. Operation starts and status of the FIXED TEMP. lamp changes from flashing to illuminated.

5. Stopping operation

Press the **RUN/STOP** key for about one second. Operation stops, the FIXED TEMP. lamp goes off and the screen switches to the initial setting screen.

When you want to correct setting errors or confirm settings


When you made a mistake during setting or reconfirm settings you made, press the **FIXED TEMP.** key again to resume setting.

When you want to change set temperature during operation, press the **FIXED TEMP.** key to enter the setting mode and change the set temperature. After change has been made, press the **ENTER** key to complete the process.

4. Operating procedures

Operating procedures (quick auto stop operation)

Procedures for quick auto stop operation

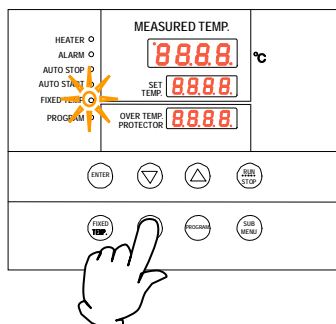

Used when you want to, for example, “stop fixed temperature operation being performed automatically in several hours. Quick auto stop operation is a function to enable auto stop timer setting during operation.

1. Setting time period before stop during fixed temperature operation

Make sure that the FIXED TEMP. lamp is illuminated to indicate the unit is the fixed temperature operation mode.

Press the **TIMER** key.

The character tim that indicates the timer is displayed on the MEASURED TEMP. display and the current set time flashes on the SET TEMP. display.

Set a time you want using the **▼▲** keys.


About the timer function

The maximum time that can be set for the timer is 999 hours 50 minutes.

Up to 99 hours 59 minutes, time can be set in minutes.

One hundred hours and over are set only in 10 minutes.

Keep the **▼▲** keys pressed to continuously change set time and you can quickly reach the time you want. Press the **▼▲** keys once at a time for fine adjustment.

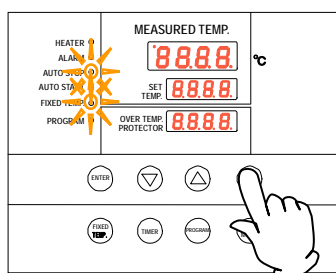

2. Starting timer operation

When you have set a time you want, start the **RUN/STOP** key for about one second. Timer operation starts with the FIXED TEMP. lamp and the AUTO STOP lamp are illuminated. Timer starts counting when the **RUN/STOP** key is pressed.

3. Stopping and ending timer operation

Operation stops automatically when the set time comes.

The buzzer sounds for about five seconds to indicate operation has stopped. At this time, the character End that indicates operation end is displayed on the SET TEMP. display with the FIXED TEMP. lamp and the AUTO STOP lamp are illuminated.

Press the **RUN/STOP** key for approx. one second to end the timer operation mode. The screen switches to the initial setting screen.

4. Operating procedures

Operating procedures (quick auto stop operation)

Change the set temp. and set time, confirm the set value

If need to change the set temp. during quick auto stop operation, press FIXED TEMP key to enter the setting mode.

If need to change the set time during quick auto stop operation, press TIMER key to enter the setting mode.

- ① After changing the time, press TIMER.
The remaining time on the timer is the time subtracting the elapsed time.
For example: set timer as 1hr30min, conduct quick auto stop operation, after 20min, change the set time to be 2hr, set it by TIMER, the remaining time is 1hr40min.
- ② After changing the time, press START/STOP for 1sec.
The quick auto stop operation proceeds again as per the changed time.
For example: set timer as 1hr30min, conduct quick auto stop operation, after 20min, change the set time to be 2hr, press START/STOP to set, the remaining time is 2hr.



Press ▼ key, switch over to display the set temp., operation mode and remaining time of set temp. display.

4. Operating procedures

Operating procedures (auto stop operation)

Procedures for auto stop operation

Used when you want to “set automatic stop after set time has elapsed from the start of the fixed temperature operation.”

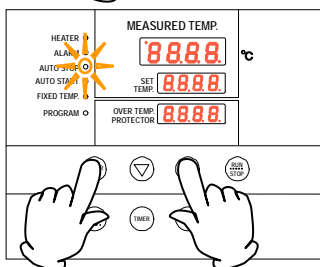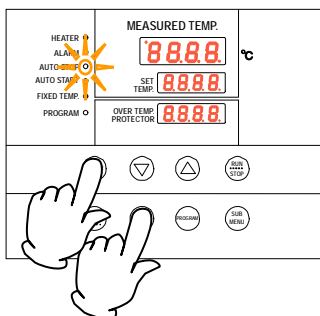

1. Setting a stop time

Press the **TIMER** key in the initial screen.

The timer mode you used in the previous session is displayed on the SET TEMP. display.

Pressing the **TIMER** key again to flash the timer mode.

Pressing the **TIMER** key again to flash the next timer mode.

Select the character AStP **AStP** that indicates the auto stop operation and press the **ENTER** key.

The character Sv **Sv** that indicate a set temperature is displayed on the MEASURED TEMP. display on which the current set temperature flashes and the AUTO STOP lamp flashes.

Set a temperature you want using the **▼▲** keys.

Press the **ENTER** key. The character tim **tim** that indicates the timer is displayed on the MEASURED TEMP. display and the current set time flashes on the SET TEMP. display.

Set a time you want using the **▼▲** keys.


About the timer function

The maximum time that can be set for the timer is 999 hours 50 minutes.

Up to 99 hours 59 minutes, time can be set in minutes.

One hundred hours and over are set only in 10 minutes.

Keep the **▼▲** keys pressed to continuously change set time and you can quickly reach the time you want. Press the **▼▲** keys once at a time for fine adjustment.

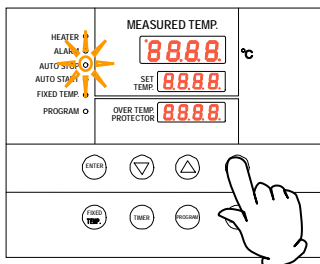

2. Starting timer operation


When you have set a time you want, start the **RUN/STOP** key for about one second.

Timer operation starts with the AUTO STOP lamp is illuminated.

Timer starts counting when the temperature in the bath (measured temperature) reaches the set temperature.


4. Operating procedures

Operating procedures (auto stop operation)

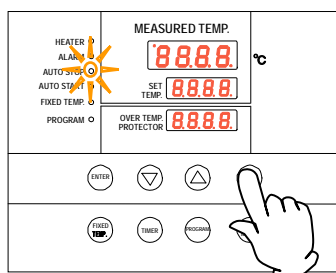

3. Stopping and ending timer operation


Operation stops automatically when the set time comes.

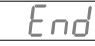The buzzer sounds for about five seconds to indicate operation has stopped. At this time, the character End  that indicates operation end is displayed on the SET TEMP. display with the AUTO STOP lamp is illuminated. Press the RUN/STOP key for approx. one second to end the timer operation mode. The screen switches to the initial setting screen.

When you want to correct set temperature or set time, or confirm settings

When you want change the set temperature or a set time, press the **TIMER** key during operation, set temperature and time for the auto stop operation with the ▼▲ keys and press the **ENTER** key to complete.

Note, however, when you change the set time you need to set a time calculated by adding the time already passed to the time to be added. Pressing the ▼ key during operation will display the set temperature, the operation mode and the remaining time on the SET TEMP. display.

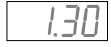
In terms of the remaining time display  a blinking dot indicates count down and an illuminating dot indicates a wait status (while temperature is increasing or decreasing to the set temperature) during which the timer has stopped counting.

4. Operating procedures

Operating procedures (auto start operation)


Procedures for auto start operation

Used when you want to “start operation automatically after the set time.”

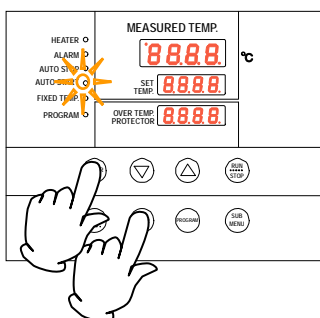

1. Setting an operation start time

- ① Press the **TIMER** key in the initial screen.
- ② The timer mode you used in the previous session is displayed on the SET TEMP. display. Pressing the **TIMER** key again to flash the timer mode. Pressing the **TIMER** key again to flash the next timer mode. Select the character AStP **AStP** that indicates the auto start operation and press the **ENTER** key.

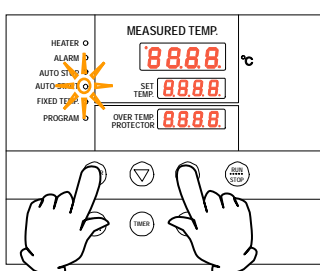

The character Sv **50** that indicate a set temperature is displayed on the MEASURED TEMP. display on which the current set temperature flashes and the AUTO START lamp flashes.

- ③ Set a temperature you want using the ▼▲ keys.
- ④ Press the **ENTER** key. The character tim **tim** that indicates the timer is displayed on the MEASURED TEMP. display and the current set time flashes on the SET TEMP. display and the AUTO START lamp flashes.
- ⑤ Set a time you want using the ▼▲ keys.


About the timer function

The maximum time that can be set for the timer is 999 hours 50 minutes.

Up to 99 hours 59 minutes, time can be set in minutes.

One hundred hours and over are set only in 10 minutes.

Keep the ▼▲ keys pressed to continuously change set time and you can quickly reach the time you want. Press the ▼▲ keys once at a time for fine adjustment.

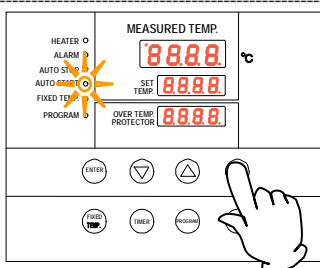


2. Starting timer operation

When you have set a time you want, start the **RUN/STOP** key for about one second.

Timer operation starts with the AUTO START lamp is illuminated.

4. Operating procedures

Operating procedures (auto start operation)

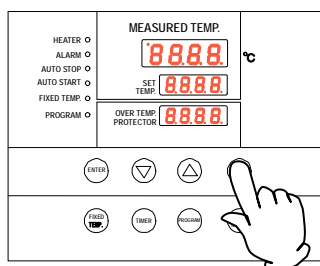

3. Stopping and ending timer operation

Operation starts automatically when the set time comes.

Press the **RUN/STOP** key for about one second to start or stop operation. The screen switches to the initial setting screen.

When you want to correct set temperature or set time, or confirm settings

When you want change the set temperature or a set time, press the **TIMER** key during operation, set temperature and time for the auto start operation with the **▼▲** keys and press the **ENTER** key to complete.

Note, however, when you change the set time you need to set a time calculated by adding the time already passed to the time to be added.

Pressing the **▼** key during operation will display the set temperature, the operation mode and the remaining time on the SET TEMP. display.


When operation has started after the auto start time, you cannot change the set time. In this case, first stop operation with the **RUN/STOP** key and repeat all settings.

4. Operating procedures

Preparing a program

Program operation

Used when you want to “operation at an increased or decreased temperature according to specific time.”

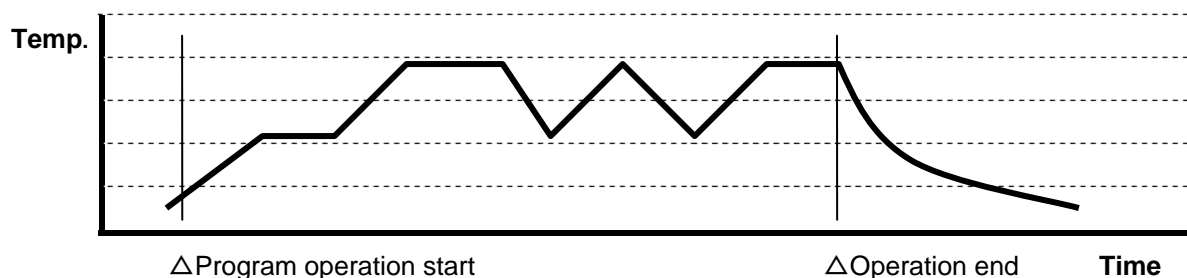

Preparing type

Up to six patterns of programs can be stored and input.

PrG1	—	One pattern of a program consisting of up to 30 steps can be made.
PrG2	PA t 1	Two patterns of a program consisting of up to 15 steps can be made.
	PA t 2	
PrG3	PA t 1	Three patterns of a program consisting of up to 10 steps can be made.
	PA t 2	
	PA t 3	

Before inputting a program

A program pattern must be registered (input) beforehand to start a program operation.

- ① Confirm the number of steps, temperatures for each step, and time on the program preparation sheet on pages 29~30 in the operation manual.
- ② Confirm the heating and cooling capacity of the device. Time needs to be set within the heating and cooling capacity of the device.
For a device having a cooling or heating capacity of 3°C for every ten minutes, for example, it requires about 35 minutes to lower or raise by 10°C from the given temperature.
- ③ Make sure that the controller has an unused pattern that meets the number of steps to be programmed.
However, when you want to use the repeat function, the number of steps to be repeated can be deducted from the number of steps for the unused pattern.

Useful function

You can use the useful repeat function when you want to repeat the same program steps. See page 28 for how to use the repeat function.

4. Operating procedures

Preparing a program

Heating • cooling time of the DVS model

Rough heating and cooling times of the DVSC model are as follows: Figures show required time for each of temperature zones. [It requires about 15 minutes to raise temperature from 100°C to 150°C.] However, be sure to conduct trial operation to determine a correct time since time required for stabilizing temperature needs to be added after the specific set temperature is reached.


Condition: Room temperature: 25°C No load Exhaust port Closed 1/3 [Unit: minute]

	DVS402C		DVS412C		DVS602C		DVS612C	
	Temp. rise time	Temp. fall time	Temp. rise time	Temp. fall time	Temp. rise time	Temp. fall time	Temp. rise time	Temp. fall time
260°C	5	-	5	-	10	-	10	-
250°C	25	10	25	10	30	10	30	10
200°C	15	20	15	20	20	20	20	20
150°C	15	30	15	30	15	30	15	30
100°C	15	40	15	40	10	45	10	45
50°C	5	100	5	100	5	75	5	75

Making a program

Programming is explained using a program pattern shown below as an example here.

1. Example of program pattern

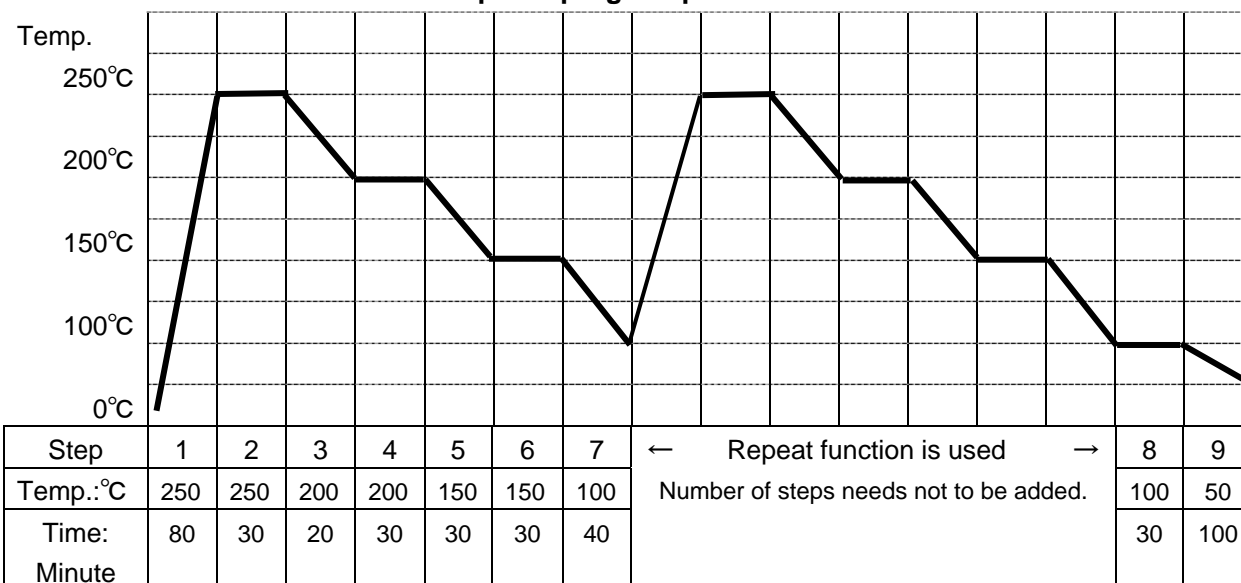

2. Turn power ON. (Turn the ELB to ON.)

When power is turned ON, the initial values will be displayed for about four seconds, then the initial screen will appear and the current bath temperature, operation mode character and the overheat prevention set temperature are displayed on each of the displays.


3. Selecting a program mode and a program pattern

- Press the **PROGRAM** key.

The program mode you used in the previous session is displayed on the SET TEMP. display.


Pressing the **PROGRAM** key again to flash the program mode.

Pressing the **PROGRAM** key again to flash the next program mode.

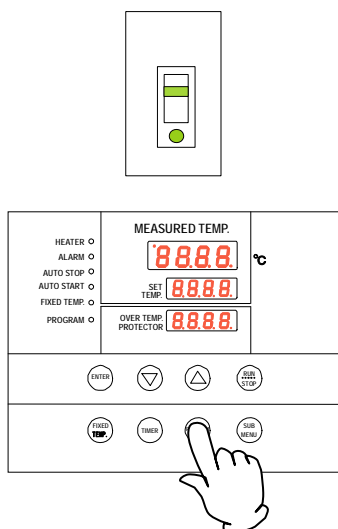

4. Operating procedures

Preparing a program

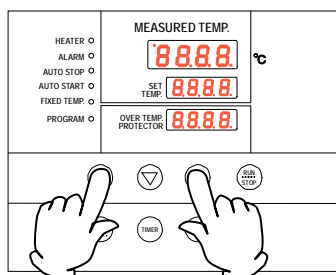


② Select a program mode you want and press the **ENTER** key.

- When PrG1 **PrG1** is selected, End **End** appears on the MEASURED TEMP. display and the number of steps registered flashes on the SET TEMP. display.
- When PrG2 **PrG2** is selected, PAT **PAT** appears on the MEASURED TEMP. display and the pattern number flashes on the SET TEMP. display. To select a pattern select [1] or [2] with the **▼▲** keys. Pressing the **ENTER** key displays End on the MEASURED TEMP. display and the number of steps registered flashes on the SET TEMP. display.
- When PrG3 **PrG3** is selected, select [1], [2], or [3] following the same procedures as for PrG2.

Since the sample program uses up to nine steps, input is allowed in any of the program modes of PrG1, PrG2, or PrG3.

Here, how to register a program is explained using PrG3 as an example.

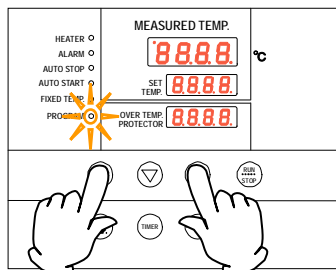

4. Registering a program (inputting a program)

① Select PrG3 following the previous section 3.

Input the number of program steps, step temperature, and step time following the entries in the programming sheet.

Press the **ENTER** key. PAT number flashes. (End will appear when PrG1 is selected. Proceed to section ⑥.)

Select an unused pattern from PAT1, PAT2, or PAT3 with the **▼▲** keys.

Press the **ENTER** key. End appears and the number of steps “n” flashes.

*End is a character that indicates the total number of steps to be used. In the example, you will input “9.”


Input the total number of program steps to be used “9” with the **▼▲** keys.


Press the **ENTER** key. The character Sv-1 **Sv-1** that indicates the set temperature of the first step is displayed and the current set temperature flashes.

Set a temperature for the first step using the **▼▲** keys.

4. Operating procedures

Preparing a program

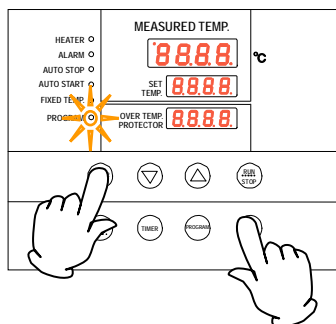

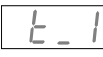Press the **ENTER** key. The character t-1  that indicates the set time of the first step is displayed and the current set time flashes.

- * To set a time you need to know the heating capacity (or cooling capacity) of the unit beforehand.
- * As an example, DVS 402C model requires about 75 minutes to increase room temperature to 250 °C . Therefore, assuming the current temperature to be 25 °C , heating capacity can be estimated roughly as three minutes to raise one °C meaning 75 minutes is necessary to increase to 250°C . In the example, the time has been estimated as 80 minutes with some allowances added.
- * The maximum time that can be set for the timer of each step is 999 hours 50 minutes.

When time has been set, press the **ENTER** key.

The temperature set character Sv-2 for the second step is displayed.

Input temperature and time of each step following the programming sheet.


When you want to repeat a certain program pattern (program repeat) in the middle of a program as in the example, a special procedure will be required. In this case, set time (t-7 in the example) for the step to be repeated (step 7 in the example) and press the **SUB MENU** key without pressing the **ENTER** key. Now you can enter the repeat function set mode.

*For operations of the program repeat function and how to register (input), follow procedures in “Program repeat operation” on page 28.

When setting the temperature and time for the final step has been completed, the screen returns to the initial setting screen.


Request for operation check

Prior to start actual operation with samples installed, be sure to perform no-load operation to check if the set temperature and the set time is appropriate.

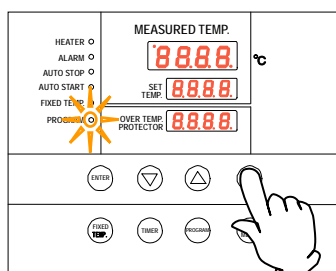

5. Starting the program operation


Press the **RUN/STOP** key for about one second. The set program operation starts.

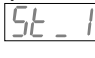The PROGRAM lamp illuminates and steps of the present operation will be displayed from the step St-1  on the SET TEMP. display.

*Press the **▼** key during operation to check the set temperature and the remaining time for each step being operated on the SET TEMP. display.

4. Operating procedures

Preparing a program

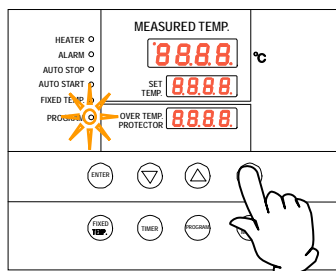

6. Terminating the program operation

When the program operation is terminated, the buzzer sounds for about five seconds to notify it to the user.

Character "End" appears on the SET TEMP. screen to indicate operation has ended.

Press the **RUN/STOP** keys to return to the initial setting screen.

About the timer function

The maximum time that can be set for the timer of each step is 999 hours 50 minutes.

Up to 99 hours 59 minutes, time can be set in minutes.

One hundred hours and over are set only in 10 minutes.

Keep the ▼▲ keys pressed to continuously change set time and you can quickly reach the time you want. Press the ▼▲ keys once at a time for fine adjustment.

When you want to correct setting errors or confirm settings

When you want to go back to previous steps for example to correct programming errors or to confirm settings, press the **FIXED TEMP.** key to return the setting screen to the previous screen.

Every time you press the **FIXED TEMP.** key you will go back by one step.

Note: Be sure to perform these operations on the program setting screen.

About wait of the program operation


When the program proceeds from a step to the next step and measured temperature has not reached or exceeded the set temperature after the set time for the step has elapsed, the next step will not start.

However, the unit has been set to proceed to the next step when the measured temperature is $\pm 3^{\circ}\text{C}$ to the set temperature.

4. Operating procedures

Program repeat operation

Using the program repeat function

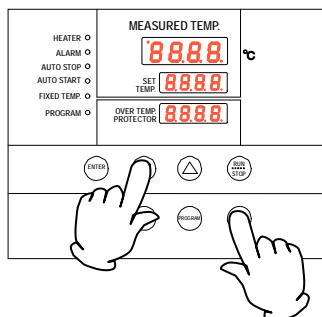

This section explains how to register a program pattern to be repeated (program repeat) during the program operation.

Using the program repeat function

This section explains how to register a program when you want to use the program repeat function in the middle of registering a program described in the previous section 4.

This procedure sets the number of step to return to “PS-n” and the number of repetition “Pc-n.” (n: step number when inputting the repeat operation)

- ① Set time (t-7 in the example) for the step to be repeated (step 7 in the example) and press the **SUB MENU** key without pressing the **ENTER** key. Now you can enter the repeat function set mode.

Character “PS-n” that indicates “return destination select” of a program pattern is displayed on the MEASURED TEMP. display. In the example, the repeat function is input in the seventh step and PS-7 **PS-7** is displayed on the MEASURED TEMP. display.

You can input a number 1 to 7 of a step to return to on the SET TEMP. display and input the number of step (1 in the example) using the **▼▲** keys.

Then press the **SUB MENU** key.

Character “Pc-n” that indicates “number of repetitions” is displayed on the MEASURED TEMP. display. Input the number (2 in the example) using the **▼▲** keys.

Pressing the **SUB MENU** key again to move the screen to the next step. The screen moves to the Sv-8 registration screen in the example.

When you want to correct setting errors or confirm settings

No alteration is allowed in the middle of the repeat set mode.

If you want to go back to the previous step for example, to correct setting errors or to reconfirm settings you have made, finish repeat setting once, press the **FIXED TEMP.** key when the screen has changed to the temperature set screen for the next step to return the set screen to the previous screen and then redo repeat setting operations.

Note: Be sure to perform these operations on the program setting screen.

If you have any questions, ask the nearest sales office or the customer support center.

4. Operating procedures

Programming sheet

Make duplicates as necessary.

Registration destination	PrG1 PrG2 PrG3 PAAt1 PAAt2 PAAt3	Control No.	
Title of test		Y/M/D	
		Prepared by	

Program pattern

															28
															27
															26
															25
															24
															23
															22
															21
															20
															19
															18
															17
															16
															15
															14
															13
															12
															11
															10
															9
															8
															7
															6
															5
															4
															3
															2
															1
700°C		600°C		500°C		400°C		300°C		200°C		100°C			分岐序号

4. Operating procedures

Programming sheet

Make duplicates as necessary.

Registration destination	PrG1 PrG2 PrG3 PAAt1 PAAt2 PAAt3	Control No.	
Title of test		Y/M/D	
		Prepared by	

Program input values



	Set temperature (°C)	Set time (unit: minutes)	Repeat function input (Return destination: No. of times)
Step 1		:	:
Step 2		:	:
Step 3		:	:
Step 4		:	:
Step 5		:	:
Step 6		:	:
Step 7		:	:
Step 8		:	:
Step 9		:	:
Step 10		:	:
Step 11		:	:
Step 12		:	:
Step 13		:	:
Step 14		:	:
Step 15		:	:
Step 16		:	:
Step 17		:	:
Step 18		:	:
Step 19		:	:
Step 20		:	:
Step 21		:	:
Step 22		:	:
Step 23		:	:
Step 24		:	:
Step 25		:	:
Step 26		:	:
Step 27		:	:
Step 28		:	:
Step 29		:	:
Step 30		:	:


4. Operating procedures

Useful functions (calibration offset function)

Using the calibration offset function

Calibration offset function compensates any differences between the target temperature in the bath and the control temperature of the controller (sensor temperature.) The function can compensate in parallel to either plus or minus side for the whole temperature band of the unit. The lock can be set or released with the **SUB MENU** keys.

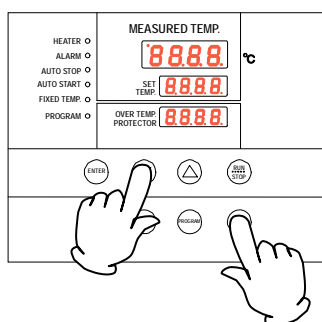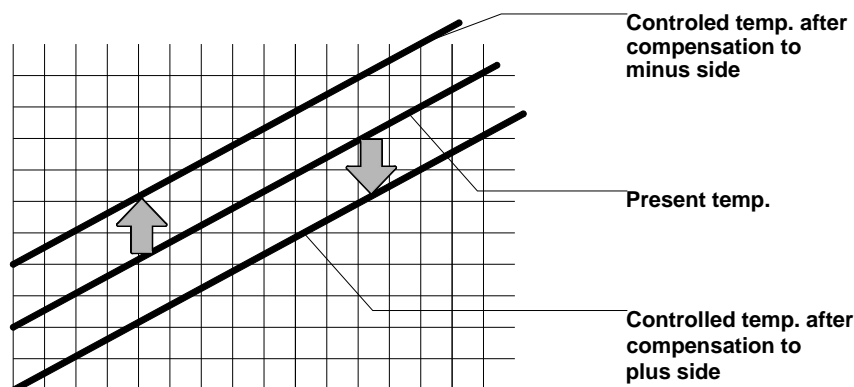


- ① Start operation at the target set temperature and confirm the temperature in the bath (sample temperature) with a temperature recorder after temperature has stabilized.
 - ② Confirm the difference between the set temperature and that in the bath (sample temperature).
 - ③ Press the **SUB MENU** key, select the character **cAL**  that indicates calibration offset using the **▼▲** keys and press the **ENTER** key.
 - ④ Input the difference between the set temperature and that in the bath using the **▼▲** keys and press the **ENTER** key to complete setting.
- * You can set either of + or – side up to +99°C and -99°C respectively for the offset compensation temperature.
When compensation is set for the – side, the MEASURED TEMP. display decreases by the compensation temperature while the temperature in the bath increases by the same amount.
When compensation is set for the + side, the MEASURED TEMP. display increases by the compensation temperature while the temperature in the bath decreases by the same amount.
 - * Since too large a compensation value may result in larger difference between the actual and indicated temperatures and may present a danger, consult our nearest sales office before entering a large compensation value.
 - * The device has, in addition to the calibration offset function, the two-point compensation function that adjusts offset for the lower temperature range and higher temperature range, for which adjustment temperatures have been input on shipping from the factory.
 - * Consult the nearest sales office before attempting validation work for the temperature adjusting device.

4. Operating procedures

Useful function (lock function)

Using the lock function

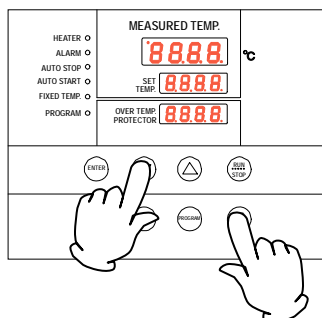

This function locks the set operation status.

The lock can be set or released with the **SUB MENU** key.

- ① Press the **SUB MENU** key, select the character Lock **LocH** that indicates setting lock using the ▼▲ keys and press the ENTER key.
- ② “OFF” is displayed on the SET TEMP. screen. To lock settings, change to “ON” using the ▲ key.
- ③ To release lock, press the **SUB MENU** key, select the character Lock **LocH** that indicates setting lock using the ▼▲ keys and press the **ENTER** key.
Lock is released when “off” is selected using the ▼ key and the ENTER key is pressed.

* When the lock function is “ON”, keys other than the **RUN/STOP** key and the **SUB MENU** key are locked.

5. Cautions on handling

Warning

1. About substances that cannot be used for the unit



Never use an explosive substance, a flammable substance or a substance containing them for this device.

An explosion or an electrical shock may result.

See section "13. List of dangerous materials" on page 44.

2. Prohibition of use/countermeasures when an error occurs



If smoke is emerges on the unit or an odd odor is felt, immediately turn the power supply off, remove the power cord (plug) from the power supply and contact your dealer or a Yamato sales office for inspection. Otherwise, a fire or an electrical shock may result. The user shall never attempt to repair the unit to avoid any possible dangers.

Caution

1. Do not place an object on the unit.



Do not place an object on the unit. Such an object may fall and cause a personal injury.

2. Do not put anything on this unit



Please do not put anything on this unit. It may cause injury. In addition, do not place inflammable materials such as paper around this unit.

3. Pay attention to high temperature scald



Due to the high temperature operation of the machine, the furnace body, the inside of the door and the samples are still in high temperature for a period of time after the operation. Pay attention not to be scalded by high temperature when taking out samples.

4. During a thunder storm



During a thunderstorm, turn off the power immediately, then turn off the circuit breaker and the main power. If this procedure is not followed, fire or electrical shock may be caused.

5. About the setting of samples



If the excessive amount of sample is set, it could be impossible to control the temperature normally. To keep the temperature control accuracy, do not use this unit in overload.

6. Return after power failure



When the operation is stopped due to power failure, after the power supply is restored, the unit can be restored to the state before the power failure. However, after the power failure, it is dangerous to run immediately. It is best to cut off the power when the power failure occurs.

5.Cautions on handling



Caution

7. Provide ventilation at the first operation



This unit exhausts smoke and smell due to burning of organic matters in the furnace when it is used for the first time. This is not abnormal, but ventilating inside the room should be done.

8. Make protective film on heater



The heater used in this unit will form a protective film on its surface under high temperature. If you usually use a relatively low temperature below 500°C, then at least rise to 700°C once, run for 10 hours to form a protective film.

9. Opening the door in high temperature will affect the unit



Please do not open or close the door when the temperature in the furnace is above 500°C. Opening or closing the door in high temperature will affect the service life of sensor, furnace body and heater. If you have to open the door, close it quickly after operation.

10. Do not leave door open for a long time at high temperature



Leaving the door open for long time at high temperature may cause a heat radiation, which could result in breakdown on the operation panel or controller.

11. Pay attention to the temperature in the furnace after the operation



After the operation, the heating stops, if the temperature in the furnace is still above 600°C, and it will take a long time to cool down to room temperature to open the door, you can turn the earth leakage circuit breaker OFF or pull out the power plug. Before opening the door, please be sure to turn on the unit again and confirm that the temperature in furnace has been lowered before opening the door to operate.

5.Cautions on handling



Caution

12. Fine powder may fly



This unit uses the iron chromium aluminum heating wire. The powder of protective film (acidized film) of heater may fly away during operation. Please protect it according to the sample condition.

13. Heater corrosion



The heater of this unit will be corroded by halogens such as chlorine and fluorine and salts contained in alkaline metals such as sodium and potassium. Do not contact the heater with such substances.

14. Sensor deterioration



The sensor used in this unit is very precise and sensitive. Please be careful not to contact with samples when taking and putting samples. Also, do not touch the sensor with bare hands, as it may deteriorate.

15. Corrosion of inner chamber and sensor stainless steel



The material of the inner chamber and sensor protection tube of this unit is SUS 310S stainless steel. It has excellent strength and oxidation resistance at high temperature, but it will be corroded by reducing substances and other reducing atmosphere at high temperature. Please do not use it.

Taking hydrogen as an example: hydrogen reduces the self-corrosion potential of SUS 310S stainless steel and promotes the anodic dissolution; hydrogen delays the formation of stainless steel passive film and decreases the corrosion resistance of passive film.

16. When using N2 gas



In the pure N2 atmosphere, the high temperature accelerates the nitriding on the surface of the heating wire and prevents the formation of the protective film. Different with atmospheric gas, it has a certain influence on its service life.

11. Notes for overheating prevention device



In case there is a small difference between the set values of temperature for overheating prevention device and that of controller, the overheating prevention device may be activated with displaying Er.19 when the temperature reaches to the set value of controller. Set the temperature of overheating prevention device to be 50°C or more higher than that of controller (when the setting temperature is lower, there is a case that overshoot occurs because this unit is high temperature drying oven. This overheating prevention device should be used to protect the unit). The default value of the overheating prevention device at factory shipment is 730°C. (Refer to P14)

6. Maintenance procedures

Daily inspection/maintenance



Warning

- Be sure to pull out the power cord for the power supply unless necessary before trying to do inspection and maintenance works.
- Start these works after the device has returned to the normal temperature.
- Never try to disassemble the unit.



Caution

- Wipe off any dirt with a tightly wrung soft cloth. Never try to clean the unit with benzene, thinner or scouring powder, or rub with a scrubbing brush. Deformation, degradation or discoloration may result.

7. When the unit is not to be used for a long time or when disposing

When the unit is not to be used for a long time or when disposing



Caution

When the unit is not going to be used for a long time

- Turn the main switch off and pull out the power cord (plug) from the power supply.



Warning

When disposing the unit

- Do not leave the unit in the area where children may have access.
- Be sure to remove handles before disposing the unit to prevent the doors from locking.
- In general, dispose the unit as a bulky waste.

Notes about disposition

Always pay attention to the preservation of the global environment.

- We highly recommend taking the unit apart as far as possible for separation or recycling to contribute to the preservation of the global environment. Major components and materials for the unit are as follows:

Names of major components	Material
Major mechanism part components	
Exterior	Steel plate SPCC (powder coating)
Furnace body	Ceramic fiber
Heat insulator	Glass wool
Door packing	Silicon rubber foam
Nameplates	Polyethylene (PET) resin film
Major electric parts	
Heater	Iron chromium aluminum electric heating wire heater
Boards	Composite parts including boards, capacitor, resistor and transformer
Power cord, wire material and others	Synthetic rubber sheathed and resin sheathed wires







8. Troubleshooting

Safety device and error codes

The unit has the self diagnostic function with a controller and a separate safety device. Table below shows possible causes and measures when the safety device is triggered.

[Error codes]

When a functional or mechanical abnormality occurs, the alarm lamp on the control panel comes on, an error code will be displayed and the alarm buzzer sounds. When an abnormality occurs, confirm the error code and immediately stop operation.

Safety device	Symptom	Possible causes and measures
Sensor error detected	Alarm lamp on  appears	<ul style="list-style-type: none"> ● Disconnection or other errors in the temperature sensor Contact our service department.
SSR short circuit detected	Alarm lamp on  appears	<ul style="list-style-type: none"> ● SSR short circuit Contact our service department.
Heater disconnection detected	Alarm lamp on  appears	<ul style="list-style-type: none"> ● Heater disconnection Contact our service department.
Memory error	Alarm lamp on  appears	<ul style="list-style-type: none"> ● Memory setting error Contact our service department.
Internal communication error	Alarm lamp on  appears	<ul style="list-style-type: none"> ● Internal communication or the temperature input circuit error Contact our service department.
Overheat	Alarm lamp on  appears	<ul style="list-style-type: none"> ● Activation of overheat prevention device Check the set temperature. Contact our service department if the unit will not recover.
Measured temperature error	Alarm lamp on ---- ---- appears	<ul style="list-style-type: none"> ● When the measured temperature is outside the displayable range Contact our service department.

8. Troubleshooting

When a malfunction is suspected

If any of the symptoms below occurs

Symptom	Check
The unit is not activated even if the power is turned on	<ul style="list-style-type: none">● Check if the power cord is connected to the power supply or to an outlet securely.● Check if power outage is occurring.
Temperature does not rise.	<ul style="list-style-type: none">● Check if the set temperature is below that in the bath.● Check if the power supply voltage has declined.● Check if the ambient temperature is low.● Check if cooling load for inside the bath is too large.
Temperature fluctuates during operation.	<ul style="list-style-type: none">● Check if the set temperature is appropriate.● Check if the power supply voltage has declined.● Check if ambient temperature fluctuates widely.● Check if cooling load for inside the bath is too large.
Displayed temperature differs from the measurement.	<ul style="list-style-type: none">● Check if the calibration offset setting is other than "0". Set it to "0." Confirm settings in "Useful functions (calibration offset function)" in page 31.

- ◆ If the symptom does not match any of the above, immediately turn the power switch on the main unit off, pull out the power cord (plug) from the power supply and contact your dealer or one of our sales offices.

9. After sales service and warranty

When requesting a repair

When requesting a repair

If any trouble occurs, immediately stop operation, turn the power switch off, pull out the power plug and contact your dealer or our sales office.

Information necessary for requesting a repair

- Model name of the product
 - Serial number
 - Date (y/m/d) of purchase
 - Description of trouble (as in detail as possible)
- } Refer to the warranty card or the nameplate on the unit.
See "3. Names and functions of parts" on page 8.

Be sure to indicate the warranty card to our service representative.

Warranty card (attached separately)

- Warranty card is given by your dealer or one of our sales offices and please fill in your dealer, date of purchase and other information and store securely.
- Warranty period is one full year from the date of purchase. Repair service for free is available according to the conditions written on the warranty card.
- For repairs after the warranty period consult your dealer or one of our sales offices. Paid repair service is available on your request when the product's functionality can be maintained by repair.

Minimum holding period of repair parts

The minimum holding period of repair parts for this product is seven years after end of production.

Repair parts here refer to parts necessary for maintaining performance of the product.

10. Specifications


Model		DR210C
Performance	Operating temperature range	300~700°C
	Temperature control precision	±5°C (at 700°C)
	Temperature rise time	Approx. 70 minutes (room temperature ~ 700°C)
Mechanism	Exterior	Cold rolled steel plate, with surface coating
	Insulating material	Aluminum silicate cotton, ceramic fiber board
	Heater	Iron chromium aluminum electric heating wire heater 1.3 kW
	Exhaust hole	I.D. 30mm x 1 on the top
Control part	Controller	VS6 program controller
	Control system	PID control of heater output with a micro computer
	Setting system	Digital setting system with the dedicated menu keys and the ▼▲ keys
	Operation modes	Fixed temperature operation, program operation, quick auto stop operation, auto stop operation, auto start operation.
	Sensor	K thermocouple
	Auxiliary functions	Lock function, calibration offset function, power outage compensation function
Safety device	Self diagnostic function	Temperature sensor error, heater error, memory error, SSR short circuit, auto overheat prevention, measured temperature error
	Protection device	Circuit Breaker, overheat prevention device (electronic type with an integrated controller)
Standard	Outer dimensions (mm) (w x d x h)	W520 × D443 × H618mm
	Inner dimensions (mm)*2 (w x d x h)	W250 × D250 × H220mm
	Internal volume	13.75L
	Weight	Approx. 24 kg
	Power supply	220V 6A
Included items		Operating instructions, warranty card

*Performance values are for the AC220V power supply with no load at an ambient temperature of 23°C.

*Operating environmental temperature range for this device is 5°C~35°C.

11. Wiring diagram

DR210C

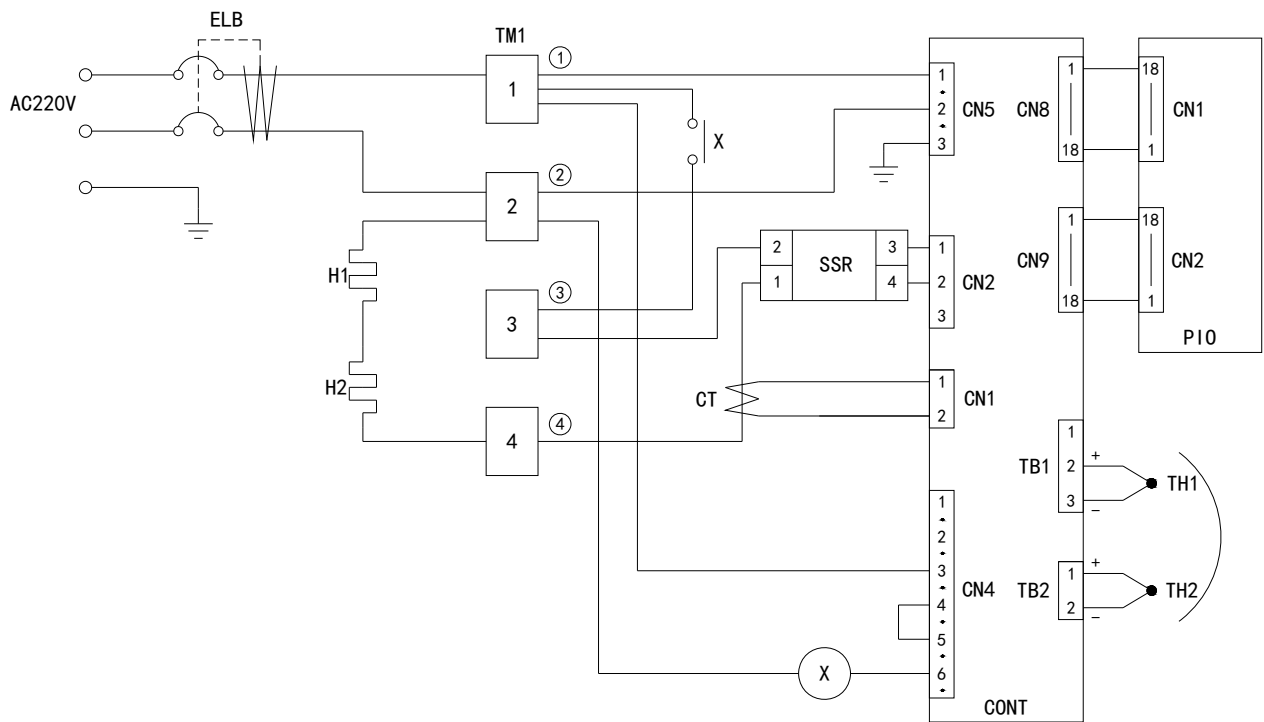

Symbol	Part name	Symbol	Part name
ELB	ELB with an over current protector	CONT	Control circuit board
TM1	Terminal block	PIO	Display circuit board
H	Heater	TH1	Control sensor
X	Main relay	TH2	Overheat prevention sensor
SSR	No-contact relay	CT	Current detection element

12. List of replacement parts

Replacement parts for DR210C

	Part name	Standard	Maker	Code No.
1	Control sensor	DR210C_01-01	Yamato	H020700004
2	VS6 Planar board	VS6	Yamato	B011401053
3	Display circuit boards	VS	Yamato	B011402007
4	Tough card	UL20861-18N*300mm	Yamato	B011299041
5	Main relay	HF116F-2/220AL1HSTFW	Yamato	A011002002
6	SSR	G3NB-225B-1 DC5~24V	Yamato	A011006011
7	Power cord kit	3*2.0sq	Yamato	A011209001
8	Circuit breaker	BV-DN IP+N 10A 30ma	Yamato	A010410007
9	Current detection element	URD CTL-6-S-4	Yamato	B010509001
10	Heater	DR210C_01-04	Yamato	H020700007

13. List of dangerous materials



Never use an explosive material, a flammable material or a material containing them for this device.

Explosive substance	Explosive substance	① Nitroglycol, glycerine trinitrate, cellulose nitrate and other explosive nitrate esters
	Explosive substance	② Trinitrobenzen, trinitrotoluenem, picric acid and other explosive nitro compounds
Flammable substances	Explosive substances	③ Acetyl hydroperoxide, methyl ethyl ketone peroxide, benzoyl peroxide and other organic peroxides
	Explosive substances	Metal "lithium", metal "potassium", metal "natrium", yellow phosphorus, phosphorus sulfide, red phosphorus, celluloids, calcium carbide (a.k.a, carbide), lime phosphide, magnesium powder, aluminum powder, metal powder other than magnesium and aluminum powder, sodium dithionous acid (a.k.a., hydrosulphite)
	Oxidizing substances	① Potassium chlorate, sodium chlorate, ammonium chlorate, and other chlorates
		② Potassium perchlorate, sodium perchlorate, ammonium perchlorate, and other perchlorates
		③ Potassium perchlorate, sodium perchlorate, ammonium perchlorate, and other inorganic perchlorates
		④ Potassium nitrate, sodium nitrate, ammonium nitrate, and other nitrates
		⑤ Sodium chlorite and other chlorites
		⑥ Calcium hypochlorite and other hypochlorites
	Flammable substances	① Ethyl ether, gasoline, acetaldehyde, propylene chloride, carbon disulfide, and other substances with ignition point at a degree 30 or more degrees below zero.
		② n-hexane, ethylene oxide, acetone, benzene, methyl ethyl ketone and other substances with ignition point between 30 degrees below zero and less than zero.
		③ Methanol, ethanol, xylene, pentyl acetate, (a.k.a.amyl acetate) and other substances with ignition point between zero and less than 30 degrees.
		④ Kerosene, light oil, terebinth oil, isopentyl alcohol(a.k.a. isoamyl alcohol), acetic acid and other substances with ignition point between 30 degrees and less than 65 degrees.
	Combustible gas	Hydrogen, acetylene, ethylene, methane, ethane, propane, butane and other Substance which is a flammable gas at 15 degrees, one air pressure.

14. Standard installation manual

*Install the product according to the following: (Confirm separately for optional items or special specifications)

Model	Serial number	Date	Installation mgr.(company name)	Installation mgr.	Judgment

No.	Item	Implementation method	TOC No. Reference page of the operating instruction manual	Judgment
Specifications				
1	Included items	Check for number of staffs against the included item field	10.Specifications field P.41	
2	Installation	• Visual check of environmental conditions Caution: Take care for environment	2. Before operating the unit • On the installation site P.4	
		•Securing a space		
Operation-related matters				
1	Source voltage	• Measure the user side voltage (outlet) with a tester • Measure voltage during operation (shall meet the specifications) Caution: Always use a plug that meets the specification for attaching to the ELB.	2. Before operating the unit • Be sure to connect the ground wire. P.4 • Power supply is P.6 10.Specifications •Specification - power supply P.41	
2	Operation start	•Starts operation Performs fixed temperature operation, auto stop operation, auto start operation, or program operation.	2. Before operating the unit P.4~7 •Installation procedures... 4. Operating procedures P.12~32	
Description				
1	Operational descriptions	Explain operations of each component according to the operational instructions	4. Operating procedures P.12~32 •Operating procedures 1. Safety precautions P.1~44 ~13.List of dangerous materials	
2	Error codes	Explain the customer about error codes and procedures for release according to the operational instructions	8. Troubleshooting ~9. After sales service and warranty P.38 40	
3	Maintenance and inspection	Explain operations of each component according to the operational instructions	6. Maintenance procedures P.36 •Daily inspection/ maintenance	
4	Completion of installation Entries	• Fill in the installation date and the installation mgr. on the nameplate of the main unit • Fill in necessary information to the warranty card and hand it over to the customer • Explanation of the route for after-sales service	9. After sales service and warranty P.40	

Responsibility

Please follow the instructions in this document when using this unit. Yamato Scientific has no responsibility for the accidents or breakdown of device if it is used with a failure to comply.

Never conduct what this document forbids. Unexpected accidents or breakdown may result in.

Note

- ◆ The contents of this document may be changed in future without notice.
- ◆ Any books with missing pages or disorderly binding may be replaced.

Instruction Manual

Hi-Temp Oven

Model DR210C

First Edition Feb. 21, 2023

Revision Jun. 28, 2023

YAMATO SCIENTIFIC CO., LTD.
Harumi Triton Square Y-36F, 1-8-11 Harumi,
Chuo-ku, Tokyo 104-6136, Japan
Tel : +81-3-5548-7122
Fax : +81-3-5548-0132
<https://www.yamato-scientific.com/>